

GETTING STARTED WITH THE CBL2 INTERFACE

The **CBL2** has a program called **DataMate** already installed. This program needs to be transferred to your calculator.

To transfer to a **TI-83 Plus** follow these instructions:

- Connect the **CBL2** to the calculator using the cable.
- Get the calculator ready to receive a program. Press **2nd**, then **LINK**. This takes you to the **SEND RECEIVE** menu. Select **RECEIVE** by using the cursor:  then **ENTER**
- Press **TRANSFER** on the **CBL2**. The **DataMate** program will now be loaded into the calculator. It takes a few minutes.
- When the transfer is complete press **2nd** and then **QUIT**
- You may now use your calculator linked to the **CBL2** interface to record and store data.

Good luck!

DataMate
successfully
installed.

Applications
menu

